

An all-Prime Editing, one-step approach for non-viral generation of a multiplex-edited CAR-T cell drug product

Emily Pomeroy, Christopher Podracky, Reyna Chang, Arika Dwivedi, Jeffrey Hussmann, Simran Padhye, Aditi Chalishazar, Kanut Laoharawee, Justin Tedeschi, Peter Chen, Alan Wilhelm, Matthew Roy, Mallikarjuna Putta, Venubabu Kotikam, Jacob Stewart-Ornstein, Seth Alexander, Joseph Elich, Hetal Patel, Andrew Anzalone, Jeremy Duffield, Jennifer Gori
Prime Medicine, Inc. 60 First Street, Cambridge, MA, USA

Background

Multiplex Prime Editing to generate an allogeneic off-the shelf cell drug product may be able to address current limitations of CAR-T cell therapy:

- Manufacturing time, costs, yield, and efficacy associated with autologous cell therapy cell quantity and quality issues
- Safety risks associated with semi-random integration & double strand breaks at multiple genomic loci

Current strategies for delivery and expression of CAR transgenes are limited by:

- Semi-random integration via lentivirus or transposons risks unintended gene disruption or activation of proto-oncogenes and may lead to variable efficacy
- Targeted integration using nuclease + template for HDR limited by low efficiency and risks associated with DSB induction (e.g., chromothripsis, p53 activation)

Limitations of current gene editing strategies for multiplex editing

- Targeted gene disruption at multiple loci simultaneously with nucleases carries a risk of chromosomal rearrangements
- Base editing to disrupt splicing or introduce pmSTOP codons is limited in scope, risks pmSTOP readthrough, and cannot support targeted integration

PASSIGE™ in combination with multiplex Prime Editing (PE) maybe be able to overcome these challenges to create a potentially best-in-class allogeneic CAR-T cell product

PASSIGE strategy to integrate CAR at T cell receptor locus (T Cell Receptor Alpha or TRAC):

- ✓ Antigen specific T cell receptors created to eliminate pathogenic cells
- ✓ Potential to prevent graft-versus-host disease by knocking out endogenous TCR expression

Prime Editing to KO MHC Class I


- ✓ Edited cells may evade patient immune system
- ✓ Allows for repeat administration if needed

Multiplex with other indication-specific Prime Edits


Prime Editing Assisted Site-Specific Integrase Gene Editing (PASSIGE):

Prime Editing in combination with integrases or recombinases for targeted integration of gene-sized DNA

- ✓ Targeted integration of DNA in a **single delivery step**
- ✓ **No double strand break (DSB)** as integrase catalyzes recombination directly
- ✓ **Integration can be irreversible:** attL and attR products are distinct from initial attB and attP sequences

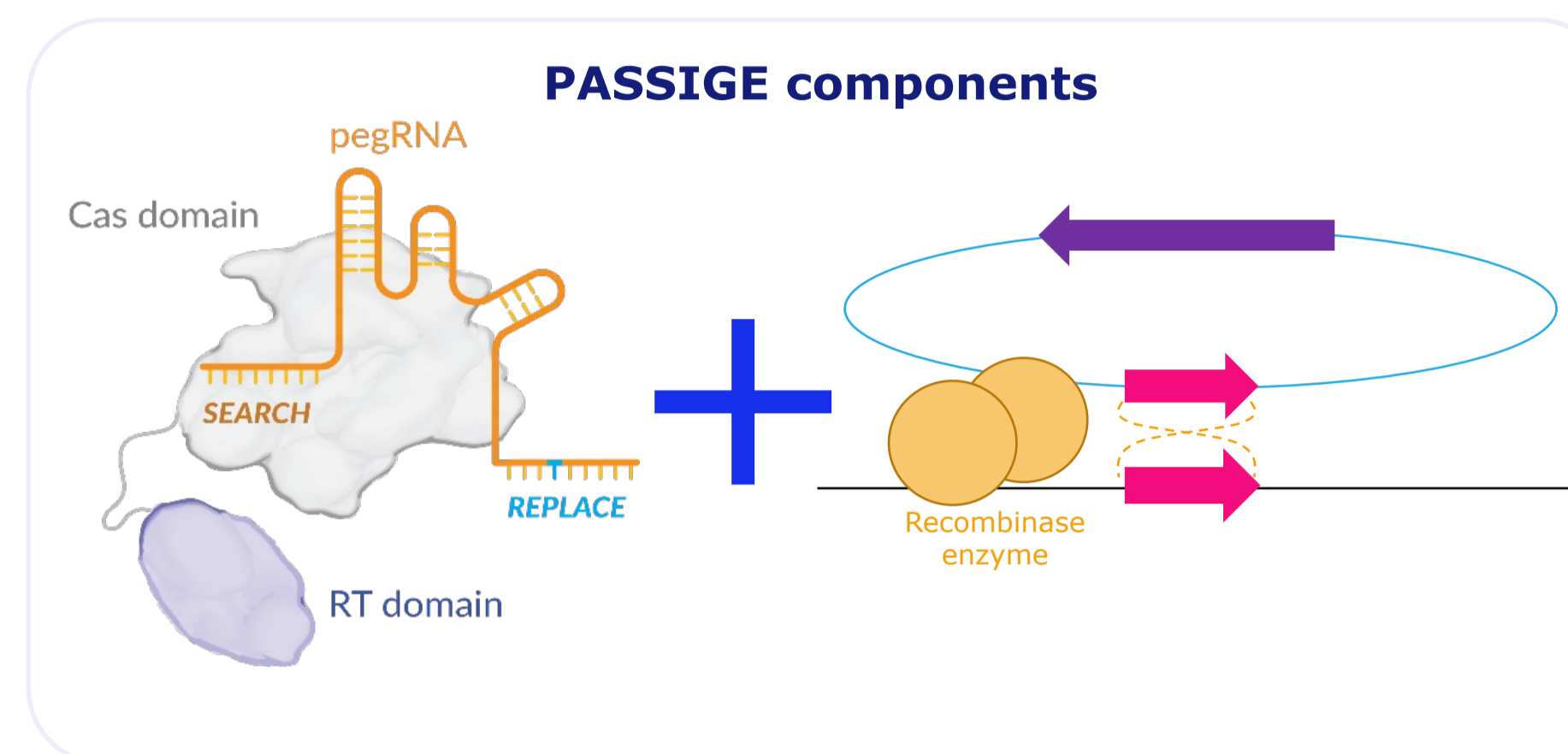

Model of Multiplex Edited CAR-T

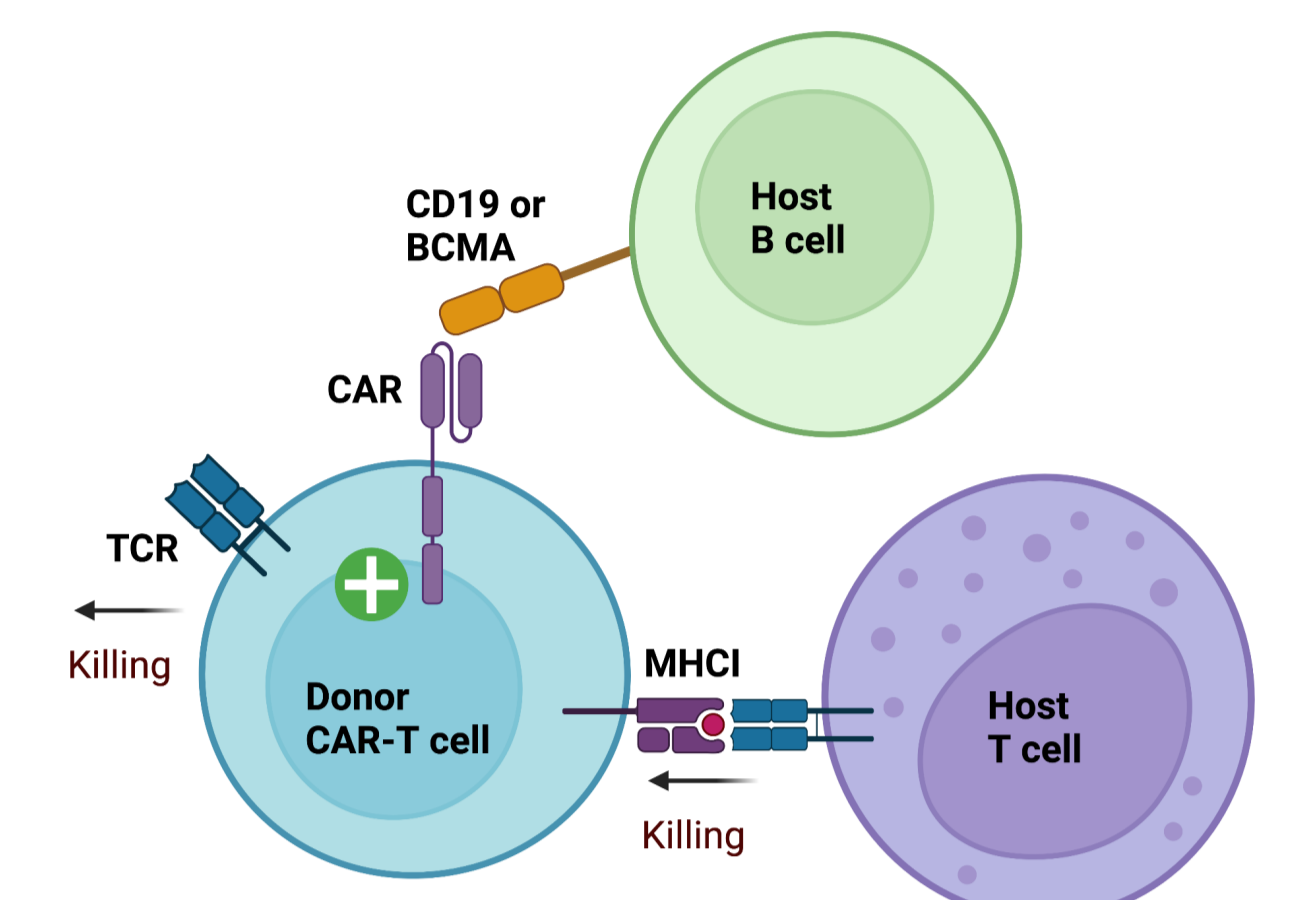

Results


1 High performing Prime Editors precisely insert attB sequence at TRAC locus with >80% efficiency

- ✓ Prime Editors screened using automated HTS in cell line
- ✓ Targeted exon 1 to make use of endogenous promoter


- ✓ Top performing Prime Editors tested in human T cells

- ✓ Top performing Prime Editors optimized in human T cells
- ✓ Achieved >80% precise attB insertion


Prime Editors screened in cell line

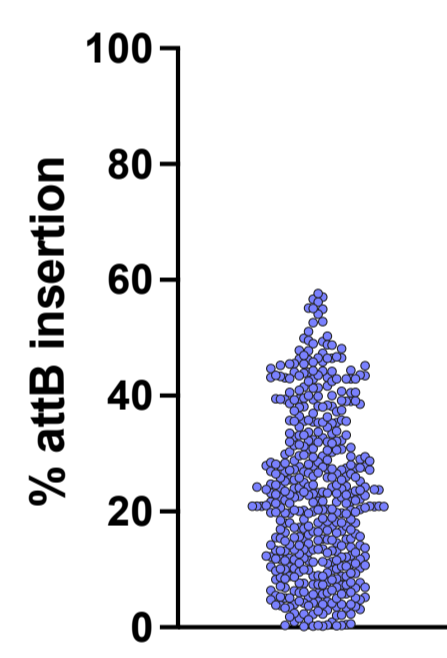

Top 3 Prime Editors validated in human T cells (n=3 donors)

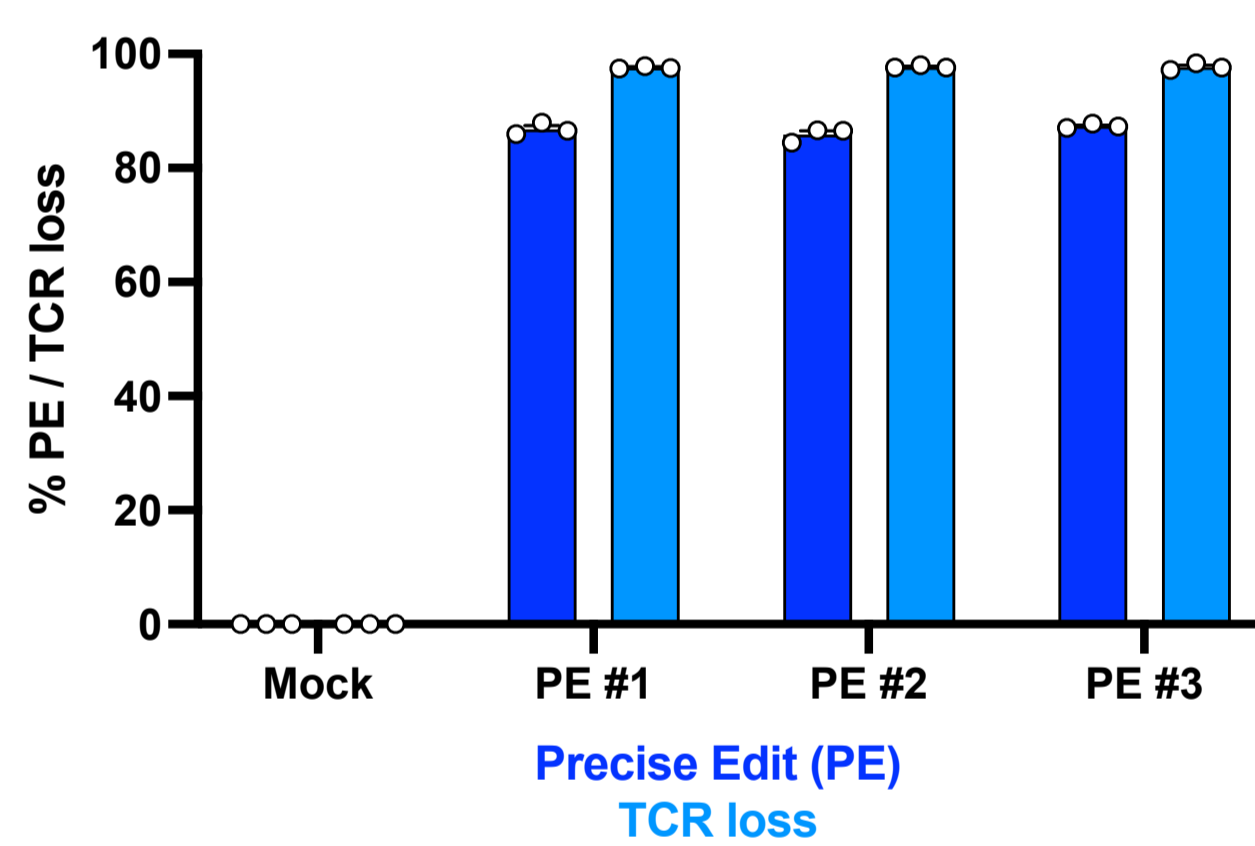


2 Non-viral PASSIGE delivery supports integration of >3.5 kb CD19 CAR construct at TRAC locus in >70% of T cells

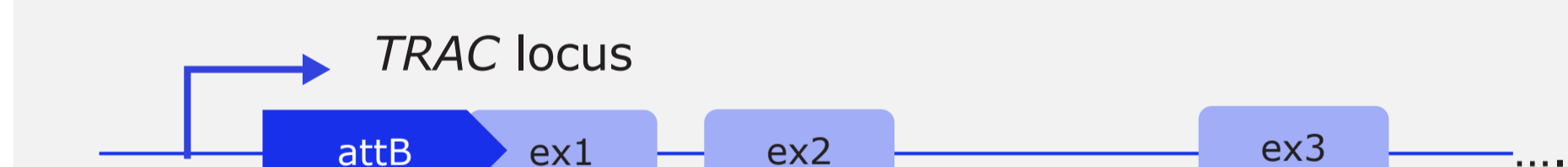

1. Prime Editing to insert recombinase sequence attB




2. Integrate CAR with Bxb1 recombinase

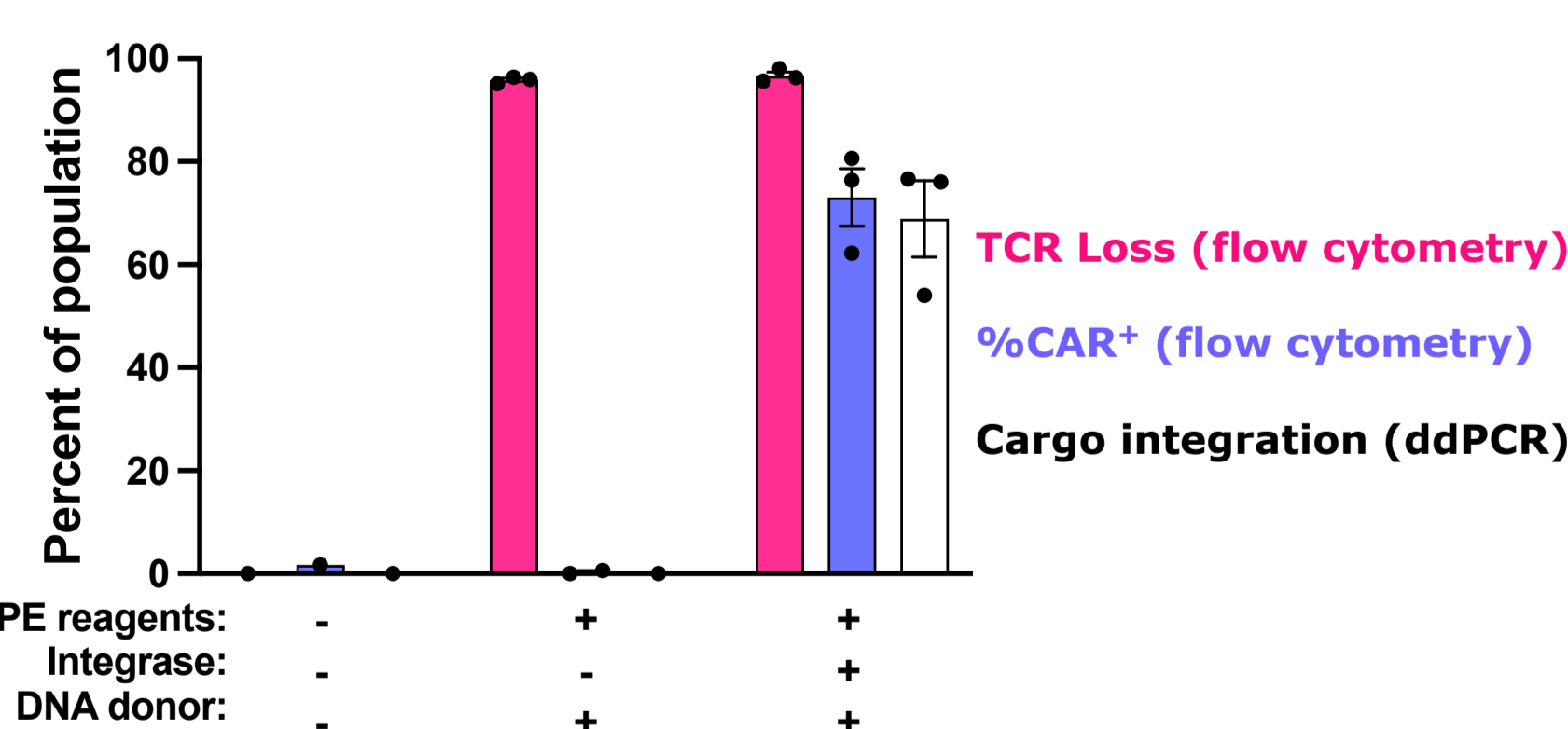

- ✓ Loss of endogenous TCR with attB insertion in TRAC exon 1
- ✓ Use of endogenous TRAC promoter allows for tuned regulation of expression
- ✓ Promoter-less cargo will not express if integrated elsewhere in genome


3 In vitro assays to evaluate CAR-T cell antigen specificity and functionality

PASSIGE CAR-T cells kill >90% of CD19+ cells

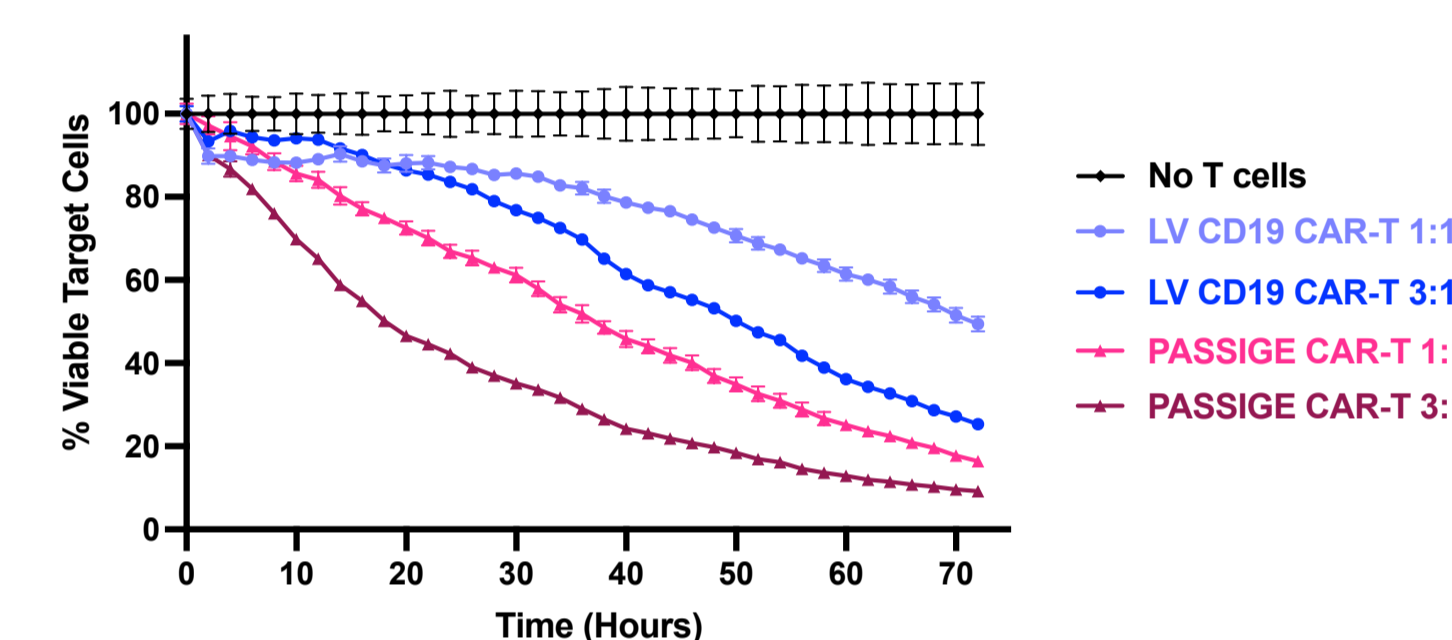


PASSIGE CAR-T: CD19-CAR at TRAC; Lentiviral vector (LV) CAR-T: CD19-CAR. PASSIGE- and LV-generated cells were ~30% CAR+ at time of plating.

PASSIGE CAR-T cells produce inflammatory cytokines in response to CD19+ cells

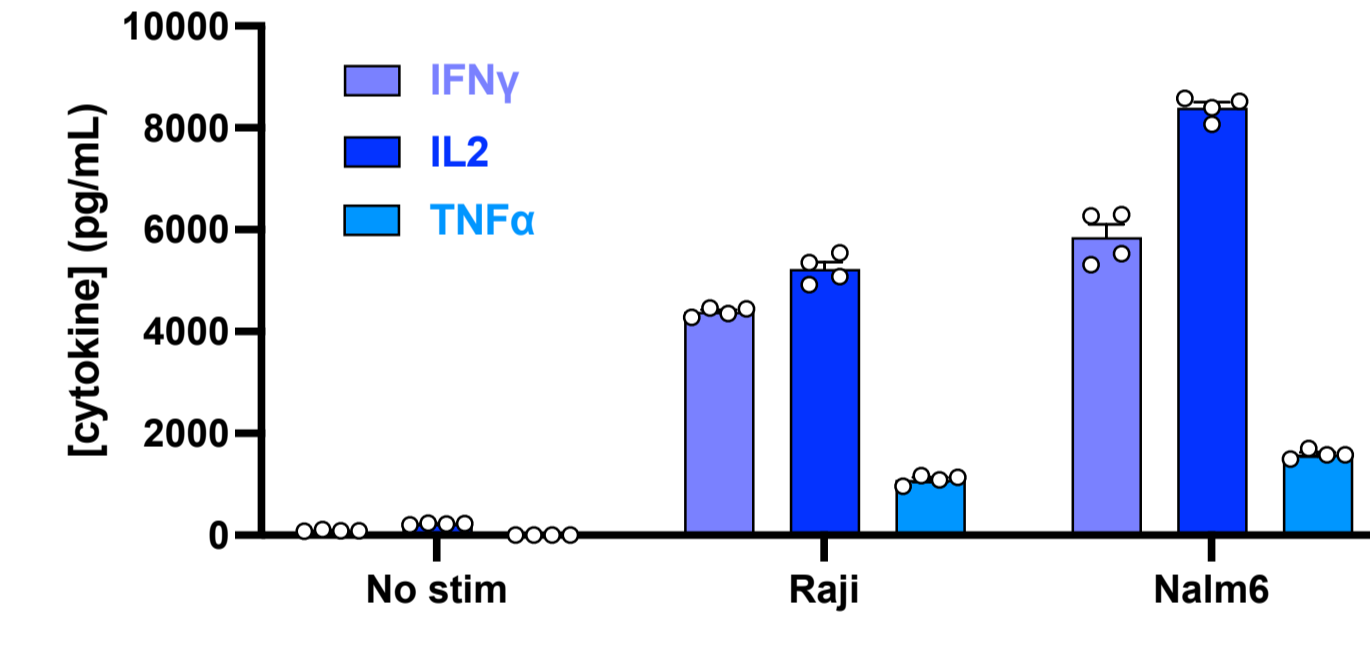


PASSIGE CAR-T: MND-CD19-CAR at TRAC; CAR-T cells co-cultured with CD19+ tumor cell lines (Raji or Nalm6) for 24 h

4 Expression of all 3 reporter cargos is consistent across DNA donor sizes (n=3 T cell donors)

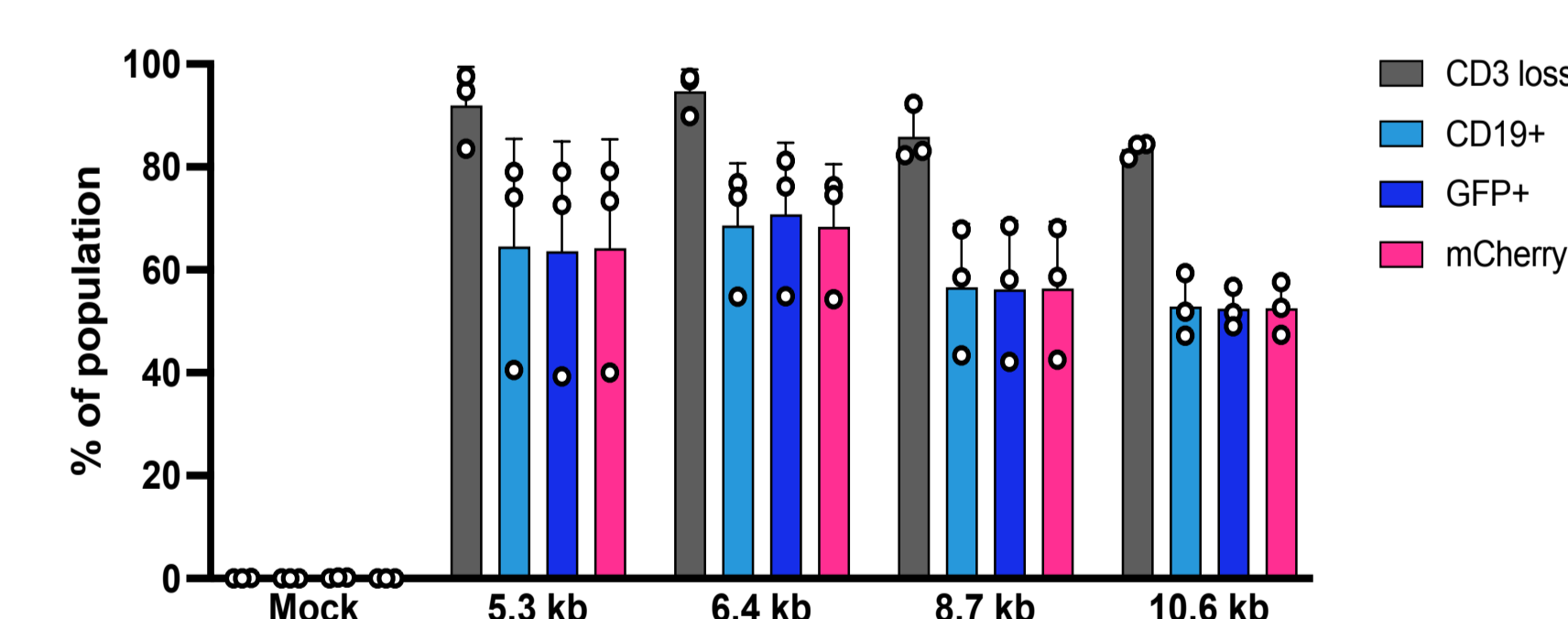

T cell health (i.e., viability and fold-expansion) is maintained regardless of DNA donor size

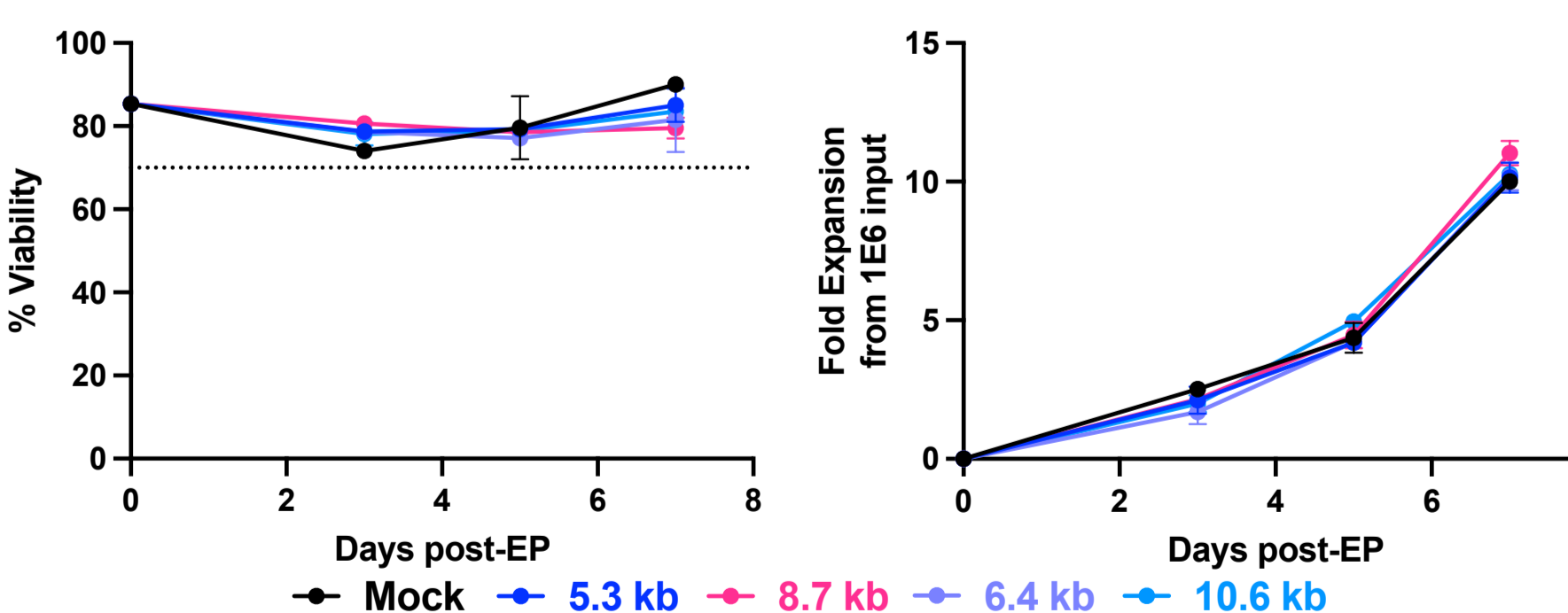


5 High performing Prime Editors precisely disrupt B2M expression with >90% efficiency in T cells

- ✓ Prime Editors screened in cell line
- ✓ Targeted early exon coding region
- ✓ Compared efficacy of multiple edit types (e.g., insertion, deletion, STOP codon)


- ✓ Top performing Prime Editors tested in human T cells

- ✓ Top performing Prime Editors optimized in human T cells
- ✓ Achieved >93% B2M protein loss

Top 3 Prime Editors validated in human T cells (n=2 donors)

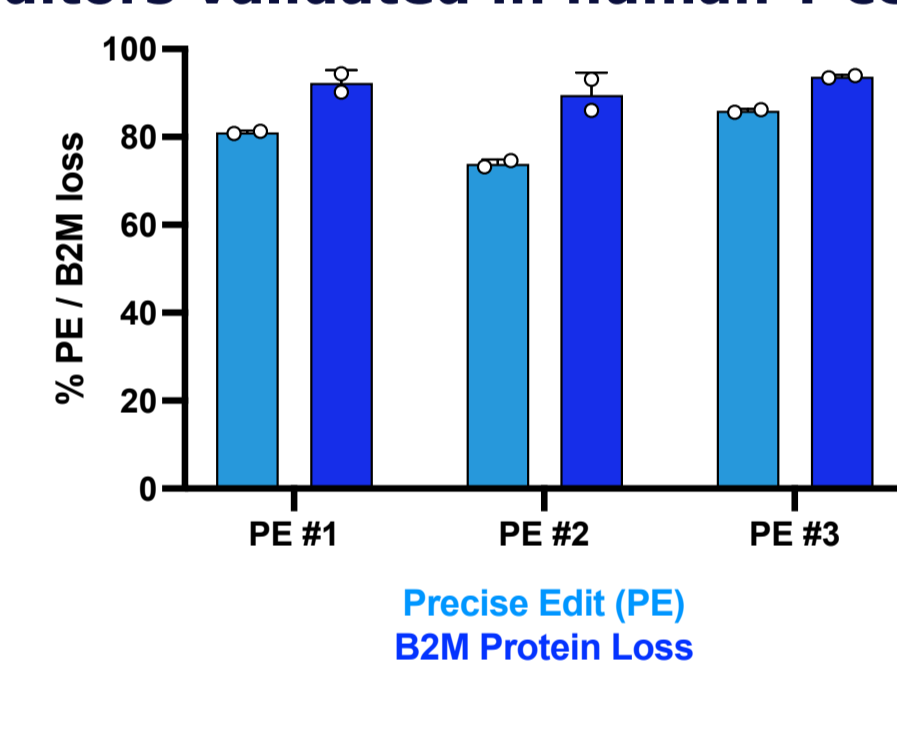


Top performing Prime Editors knock out B2M protein expression in >90% of T cells

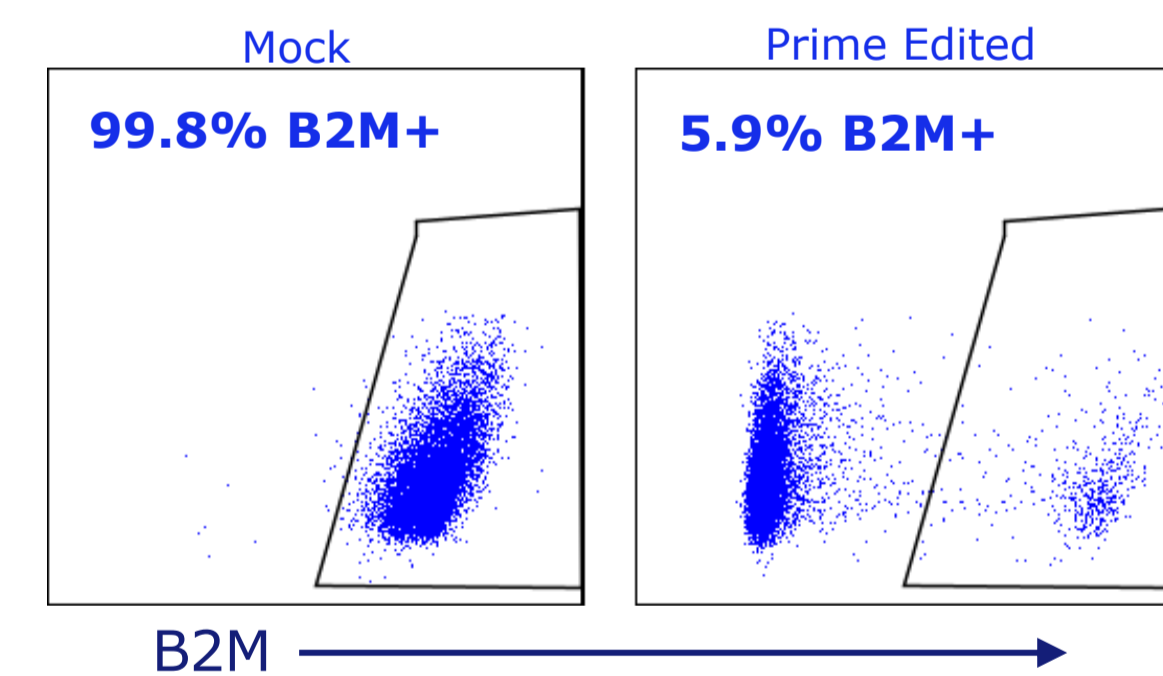

6 Prime Editing efficiency is maintained in multiplex context at 5 target sites in T cells

Percent precise Prime Editing in multiplex normalized to Prime Editing efficiency in singleplex for each target site.

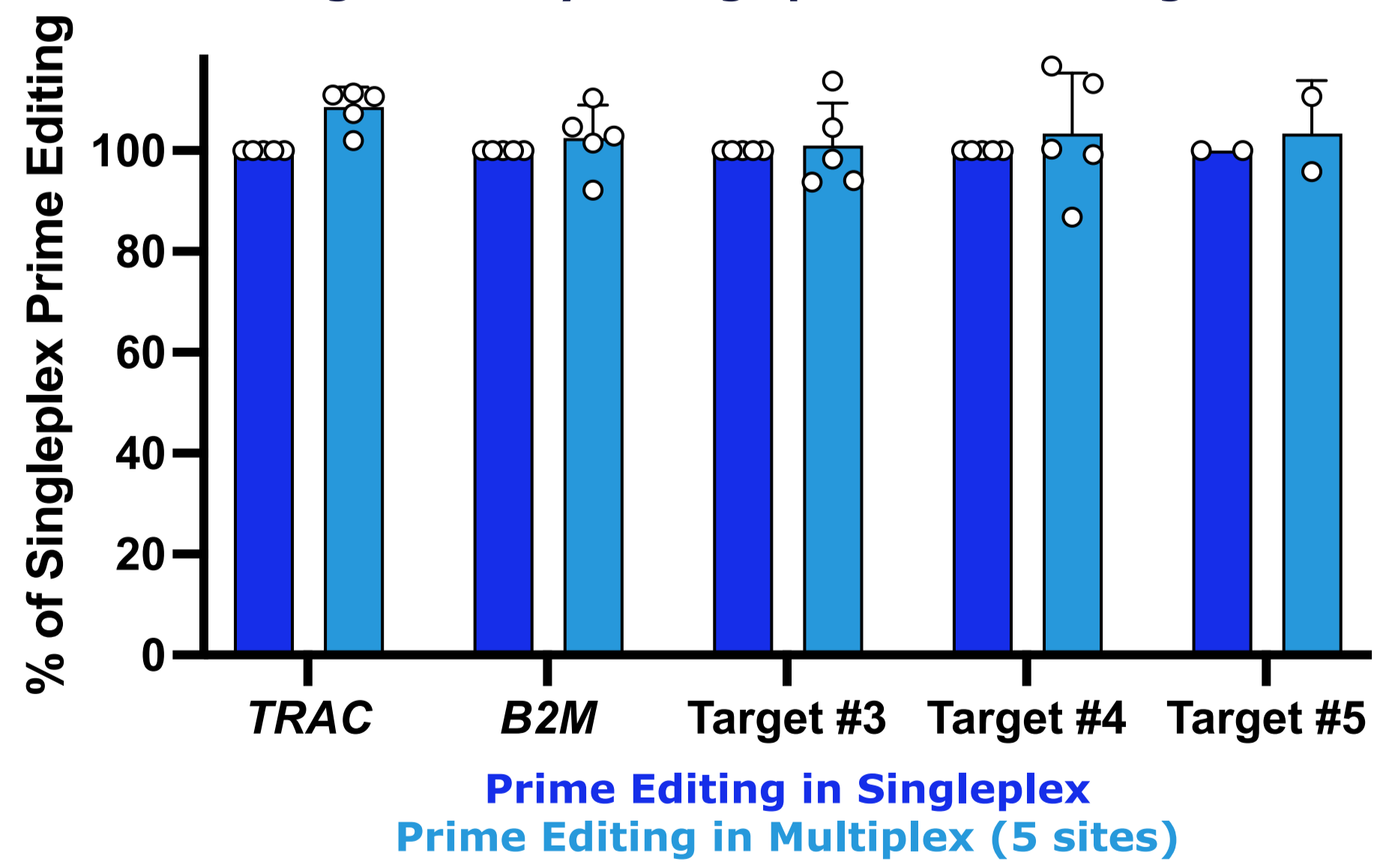

Conclusions

- Prime Editing precisely introduced recombinase target sequence at TRAC locus in >80% of T cells
- Achieved >70% site-specific integration of CD19-targeting CAR through systematic PASSIGE component and process optimization
- PASSIGE can be multiplexed with Prime Editing at other target sites in a single delivery step with no loss of efficiency observed
- PASSIGE-generated CAR-T cells are healthy and show potent antigen-specific function and cytotoxicity